


Rotary Valves

MC6000 Replacement Valves - 1/4-28 Thread Ports

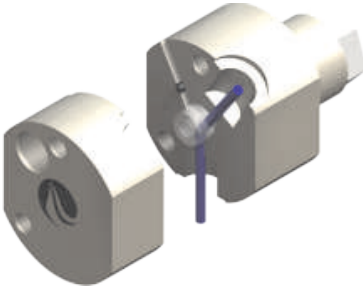

Note: Valve body shown sectioned for clarity.

Description	Materials (Body/Plug)	Catalog Number	Flow Path (Blue)
"Y" Valve 120° 3-port	PEEK/UHMWPE	10153	

VALVE REPLACEMENT

To replace the valve, prime all fluid from the syringe, and follow the steps below:

1. Remove syringe.
2. Disconnect the tubing from the valve.
3. Remove the two valve screws used to secure the valve to the panel.
4. Remove the valve from the pump.

To install the valve follow the below steps:

5. Align internal valve "D" slot with the flat on the motor shaft.
6. Gently push and rotate the valve into the correct position to align the pins with the holes in the panel.
7. Securely tighten (clockwise) the plunger lock screws.
8. Initialize the pump.

